



WATERBORNE™, INC.

Clinical & Environmental Parasitology Products ♦ 6045 Hurst Street, New Orleans, LA, USA 70118-6129
Tel. & FAX: 504-895-3338 or Toll-Free: 866-895-3338 ♦ Email: custserv@waterborneinc.com ♦ Website: www.waterborneinc.com

M101. No-Fade™ Mounting Medium

This product consists of 3.5 mL of mounting medium containing the anti-fading agent, DABCO, in glycerol with formalin. No-Fade™ Mounting Medium has been created for use with fluorescein-labeled antibodies and is conveniently packaged in a plastic dropper vial. This product can be used to mount 70-75 slide wells, using one drop (~ 45 uL) per well. The shelf-life of this product is approximately 1 year, when stored at 4° C.

NOTE: Store at 4° C and minimize exposure to light. Some yellowing may occur over time with exposure to light - this will not affect performance.

M102. Elvanol No-Fade™ Mounting Medium

This product consists of 3.5 mL of mounting medium consisting of polyvinyl alcohol, DABCO (anti-quenching agent), and glycerol in water. It is conveniently packaged in a plastic dropper vial and can be used to mount approx. 75 slide wells, using one drop per well. After mounting a slide with this medium, one may view the slide immediately and/or allow the medium to dry overnight at 25° C or 37° C to form a firmer, more durable preparation for long-term storage and repeated viewing. After drying, keep slides in the refrigerator or the freezer in a sealed, light-proof container for long-term storage.

NOTE: Store at room temperature wrapped in aluminum foil to protect from light. If, during storage, the medium forms a gel, simply heat the vial in a beaker of hot water for several minutes, or microwave with the cap loose for 7-10 seconds, to restore it to a liquid.